



BeadLogic installation guide

Windows


Download and double click the application (.exe) file from the Internet (or double click the .exe file in the dongle) to begin the installation.

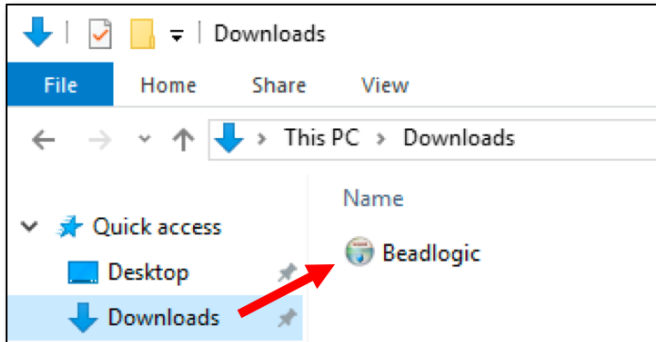

Follow the steps in the BeadLogic Setup wizard.

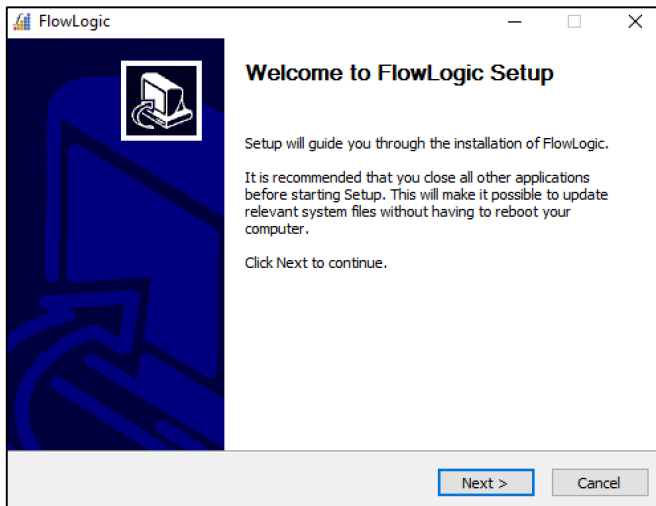

Launch BeadLogic from the applications menu or from the shortcut on the desktop. BeadLogic can also be uninstalled using the option in the applications menu.

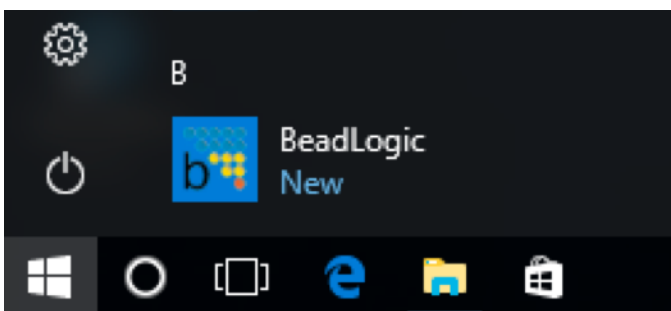

Apple

Download and double click the BeadLogic disk image (.dmg) file from the Internet (or double click the .dmg file in the dongle) and drag the BeadLogic application into the Applications folder.



When launching BeadLogic for the first time, a message may appear stating that BeadLogic can't be opened because it is from an unidentified developer. In order to run the program, right click on the application icon and then choose 'Open', and again 'Open' from the subsequent window.



Once BeadLogic has been run once, simply double click the Beadlogic application icon in the Applications folder or in Launchpad to launch the program.



Running BeadLogic from the dongle

For optimal speed and performance, it is recommended to install BeadLogic to the computer. However, it is possible to run directly from the dongle.

To run from the dongle, open the folder titled 'BeadLogic run from USB'. For Windows and Linux users, double click the .jar file to launch the program. For Apple users, double click the BeadLogic application file.

Please be aware that a more recent BeadLogic version update may exist. If you wish to run BeadLogic offline but with the current software version, take the dongle to a computer that is online, download the latest version from the Miltenyi Biotec website and replace the original application on the dongle with the new download. This can now be run directly from the dongle or copied and pasted to replace the version on other computers.

As new features are added with software updates, it may be the case that analysis files saved with newer versions of BeadLogic may not open correctly with older versions. Please also note that the licence contained in the dongle is linked to the hardware and so will not work if transferred to a new location.

